



## Bio Pump Operation Guide

- 1) Turn on pump utilizing the Power button.



- 2) Press Quick Sample until "5 Minutes" appears on the screen.



- 3) Place the Air-O-Cell cassette on the inlet of the pump, with the arrow pointing down.



- 4) Press Start/Stop to begin sampling.



- 5) When sampling is complete, the Bio Pump will beep and flash a yellow light. You may remove your cassette.



For more instructions on handling the Air-O-Cell cassette, please reference the Microbial Sampling Procedures that will be located in your paperwork from this order.