



QuickTake 15 Pump Instructions

1. Place pump in desired location.
2. Place Air-O-Cell cassette on the rubber grommet on top of pump so it fits snugly and air flow direction arrow points towards pump.
3. Simultaneously press and hold up and down arrow at same time until LED screen will lights up and will display “t 5”. The number indicates the pump run time.
4. Simultaneously press and release up and down arrow and pump will start running.
5. LCD screen should show “t _” where the _ will be the number of minutes remaining in sample time.
6. Once sampling is completed, the unit will beep 3 times, and display will show “done”.
7. If you are doing multiple samples, simultaneously press up and down arrows, “t 5” will show on LED. Replace Air-O-Cell cassette, and press and release up and down arrows again to start pump. Repeat until all cassettes have been sampled.
8. Remove cassette from pump by gently pulling upward. Pump will automatically shut off after 5 – 10 minutes of non-use.
9. Seal cassette, document sample run time and return to SGS Galson for analysis.

Charging the equipment

1. Plug charger into port on side of pump.
2. Charge indicator light will illuminate on bottom right of LED screen.
3. A full battery charge takes approximately 5 hours.

If you have any questions regarding this pump or the sampling process, please feel free to contact SGS Galson at (888) 432-5227

GALSON LABORATORIES INC. IS NOW PART OF SGS, THE WORLD'S LEADING INSPECTION, VERIFICATION, TESTING AND CERTIFICATION COMPANY.