



GALSON

Removing Custody Seal for Sampling

- Place finger on cassette nib, use provided screwdriver to perforate the seal in circular motion around the inside of the cassette's top ring.



- Remove the perforated piece, discard, and repeat with the bottom of the cassette.



- If sampling open-faced or using aluminum cyclone, perforate the seal in the first groove where the top ring meets the middle ring. Remove the perforated piece, discard, and repeat with the bottom of the cassette.



NOTICE: Be sure to clear your nibs of all debris prior to reinserting into cassettes. Failure to do so may cause debris to enter your sample.

GALSON LABORATORIES INC. IS NOW PART OF SGS, THE WORLD'S LEADING INSPECTION, VERIFICATION, TESTING AND CERTIFICATION COMPANY.